



Tables in JASON

a Cheat Sheet




Tables in JASON come in many forms including Parameter, Peak, Integral/Multiplet and Arrayed tables.

The properties of a table can be changed in the Table Tools Panel.

Properties which can be changed include:

- Setting the currently selected options as the default for all new tables created.
- Displaying the row and column indexes.
- Selecting the font for the table column headers.

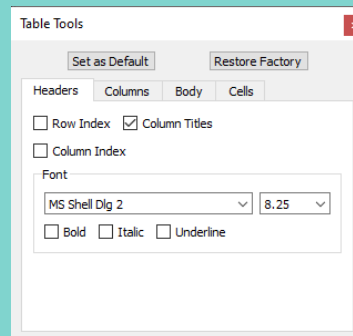

	Columns	Units	Decimal Places	Alignment	Filter
A	<input checked="" type="checkbox"/> Pos (ppm)	ppm	3	Right	n/a
B	<input checked="" type="checkbox"/> α -Pos (ppm)	ppm	5	Right	n/a
C	<input checked="" type="checkbox"/> Height	n/a	2	Right	n/a
D	<input checked="" type="checkbox"/> α -Height	n/a	4	Right	n/a
E	<input checked="" type="checkbox"/> Width (Hz)	Hz	2	Right	n/a
F	<input checked="" type="checkbox"/> α -Width (Hz)	Hz	4	Right	n/a
G	<input checked="" type="checkbox"/> Kurtosis	n/a	2	Right	n/a
H	<input checked="" type="checkbox"/> α -Kurtosis	n/a	4	Right	n/a
I	<input checked="" type="checkbox"/> Area	n/a	2	Right	n/a
J	<input checked="" type="checkbox"/> α -Area	n/a	4	Right	n/a
K	<input checked="" type="checkbox"/> Type	n/a	n/a	Right	compound contaminant NMR solvent reaction solvent 13C satellite sideband
L	<input checked="" type="checkbox"/> Label	n/a	n/a	Right	

Customize table columns

- Show or hide columns using the check box.
- Change the units displayed by double clicking on the unit and selecting from the drop down list of available units.
- Set the number of decimal places
- Define the alignment in the columns
- Where available use the filter function to show only the values you want to see.
- Rearrange table columns by drag & drop
- Sort column contents by clicking on the header

Define the table body settings

- Define the font for the main table body
- Show rows in alternating colors.
- Set the font for individual cells (*This won't be saved as default*)

